


Zebra & Quagga Mussel Veliger Detection Services

VELIGER SAMPLE SHIPPING PROTOCOL (INSTRUCTIONS)

“WHAT'S IN YOUR WATER?®”

Protocol for shipping samples collected with Clean Lakes, Inc. Veliger Sample Collection Kit:

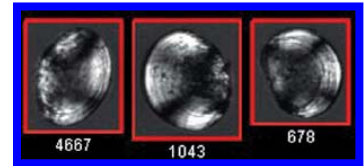

- 1: The plastic containers (bottles) provided include a premeasured amount of ethanol to preserve the sample. Fill the sample container to the full mark. **Do not overfill.**
2. Samples must be in the plastic containers (bottles) provided with screw lid securely fastened with tape. Up to **two (2)** of the sample bottles provided in the **Veliger Sample Collection Kits** can be shipped at a time.
3. Place all containers into provided sealable plastic bag (e.g. Zip Lock) and then place this into the other sealable plastic bag provided (e.g. Zip Lock).
4. Place sealed bags and sample container into the primary shipping box provided. Seal this box with clear packing tape (tape not provided).
5. Place primary shipping box into the secondary shipping box provided. Seal this box with clear packing tape (tape not provided).
6. Include a complete return address on the package, affix proper postage (or ship via Federal Express or United Parcel Service) and ship to

Clean Lakes, Inc.
Attn: Veliger Detection Division
1770 West State Street #125
Boise, ID 83702

Note: If you purchased one of CLI's **Veliger Sample Collection Kits**, sample containers and the supplies listed above for sample shipping are included. Simply take the sample bottle out of its individual packaging, collect the sample, add the sample to the sample container, and place the sample container back into the shipping package in the same way it was received. If you did not purchase one of CLI's **Veliger Sample Collection Kits**, follow the shipping protocol included with the Sample Collection Protocol.

For more information, contact Leif Elgethun, E.E.I., LEED AP, CLI's Zebra & Quagga Invasive Mussel Veliger Detection Manager, at Clean Lakes, Inc. Attn: Veliger Detection Division, 1770 West State Street #125 Boise, ID 83702 via (cell: 208-301-2293), (Fax: 888-330-8602), or (Email: lelgethun@cleanlake.com), or contact one of our offices below.

Pacific Northwest: Coeur D'Alene, Idaho: Phone: 208-665-1475, Fax: 208-665-1479
San Francisco Bay Area: Martinez, California: Phone: 925-957-1905, Fax: 925-957-1906
Santa Monica Bay Area: Westlake Village, California: Phone: 818-889-8691, Fax: 818-889-8693